

**ATCC medium: 1948 LB medium (ATCC medium 1065) with 50 mcg/ml ampicillin and 20 mcg/ml kanamycin**

ATCC Medium 1065 (see below) with 50.0 mcg/ml ampicillin and 20 mcg/ml kanamycin

Filter-sterilize solutions of ampicillin and kanamycin and add to sterile, cooled medium to achieve a final concentration of 50.0 mcg/ml and 20.0 mcg/ml respectively.

*ATCC Medium 1065:*

|                                |        |
|--------------------------------|--------|
| Tryptone (BD 211705).....      | 10.0 g |
| Yeast Extract (BD 212750)..... | 5.0 g  |
| NaCl.....                      | 10.0 g |
| Distilled water.....           | 1.0 L  |

Autoclave at 121C for 15 minutes.