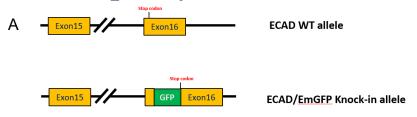


Technical Data Sheet: PANC-1 ECAD-EmGFP MET Reporter Cell Line

ATCC® Number	CRL-1469MET™
Organism	Homo sapiens, human
Tissue/Disease Source	Pancreas, pancreatic carcinoma
Product Description	The PANC-1 E-cadherin (ECAD)-EmGFP reporter cell line was created by knocking-in an EmGFP reporter gene into the endogenous E-cadherin gene using CRISPR/Cas9 technology. We confirmed that EmGFP activities accurately report E-cadherin gene expression, and showed that ECAD-EmGFP expression can be induced by microRNA-200 treatment, indicating cells transition to epithelial status.
Application	This cell line is not only a useful in vitro model for dissecting the molecular switches underlying EMT and MET, but could also be used for screening compounds targeting EMT or MET in pancreatic cancer.

E-cadherin gene expression in PANC-1 ECAD-EmGFP cells



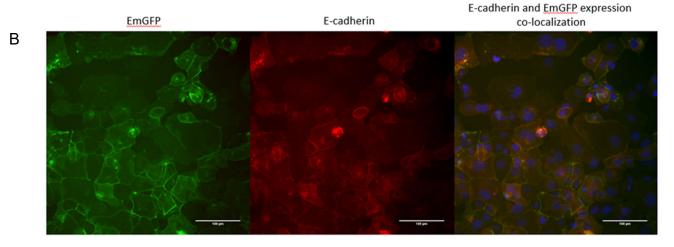


Figure 1. EmGFP activities accurately report E-cadherin gene expression in PANC-1 ECAD-EmGFP cells (ATCC® CRL-1469METTM). (A) Partial schematic diagram of the E-cadherin wild type allele and ECAD-EmGFP knock-in allele, in which EmGFP is incorporated adjacent to endogenous E-cadherin in Exon 16. (B) Endogenous EmGFP (left, green) co-localized with E-cadherin detected by immunofluorescence assay (middle, red) as shown in the merged image (right).

Cell Morphology

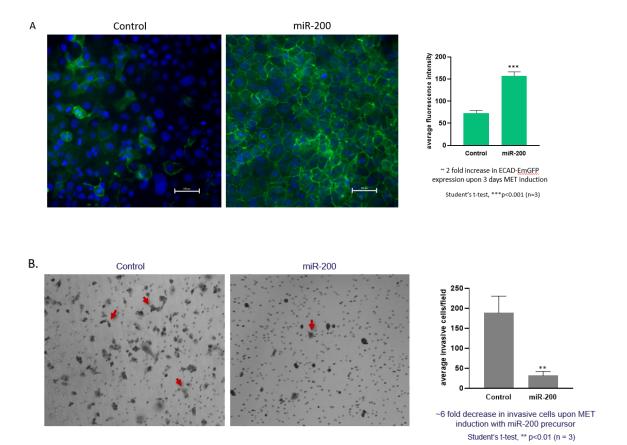


Figure 2. MicroRNA-200 treatment results in the induction of ECAD-EmGFP expression, decrease in SNAIL expression, and decrease in the invasive capabilities of PANC-1 ECAD-EmGFP cells. (A) PANC-1 ECAD-EmGFP cells were treated with miRNA-200 or an equivalent volume of 1x PBS (as a control) for 3 days. miRNA-200 treatment induced a significant increase in ECAD-EmGFP expression. The nuclei of the cells were counterstained with DAPI. (B) After 3 days miR-200 treatment, PANC-1 ECAD-EmGFP cells were monitored over a 48 hours period for invasion through a trans-well invasion assay. The induced cells showed a significant decrease in invasive capacity.

© 2020 American Type Culture Collection. The ATCC trademark and trade name, and any other trademarks listed in this publication are owned by the American Type Culture Collection unless indicated otherwise.

These products are for laboratory use only. Not for human or diagnostic use. ATCC products may not be resold, modified for resale, used to provide commercial services, or to manufacture commercial products without prior ATCC written approval.