

Streamline Your Microbial QC Testing with Confidence, Ease, and Consistency

Nilay Chakraborty, PhD, MBA
BioNexus Foundation Principal Scientist
and Head of Cryobiology, ATCC

ASM Microbe 2025
Industry & Science Showcase




About ATCC


- Founded in 1925, ATCC® is a non-profit organization with HQ in Manassas, VA, and an R&D and Services center in Gaithersburg, MD
- World's premier biological materials resource and standards development organization
- 5,000 cell lines
- 80,000 microorganisms
- Genomic & synthetic nucleic acids
- Media/reagents
- ATCC® collaborates with and supports the scientific community with industry-standard biological products and innovative solutions
- Growing portfolio of products and services
- Sales and distribution in 150 countries, 20 international distributors
- Talented team of 600+ employees, over one-third with advanced degrees




Enabling scientific progress for nearly 100 years

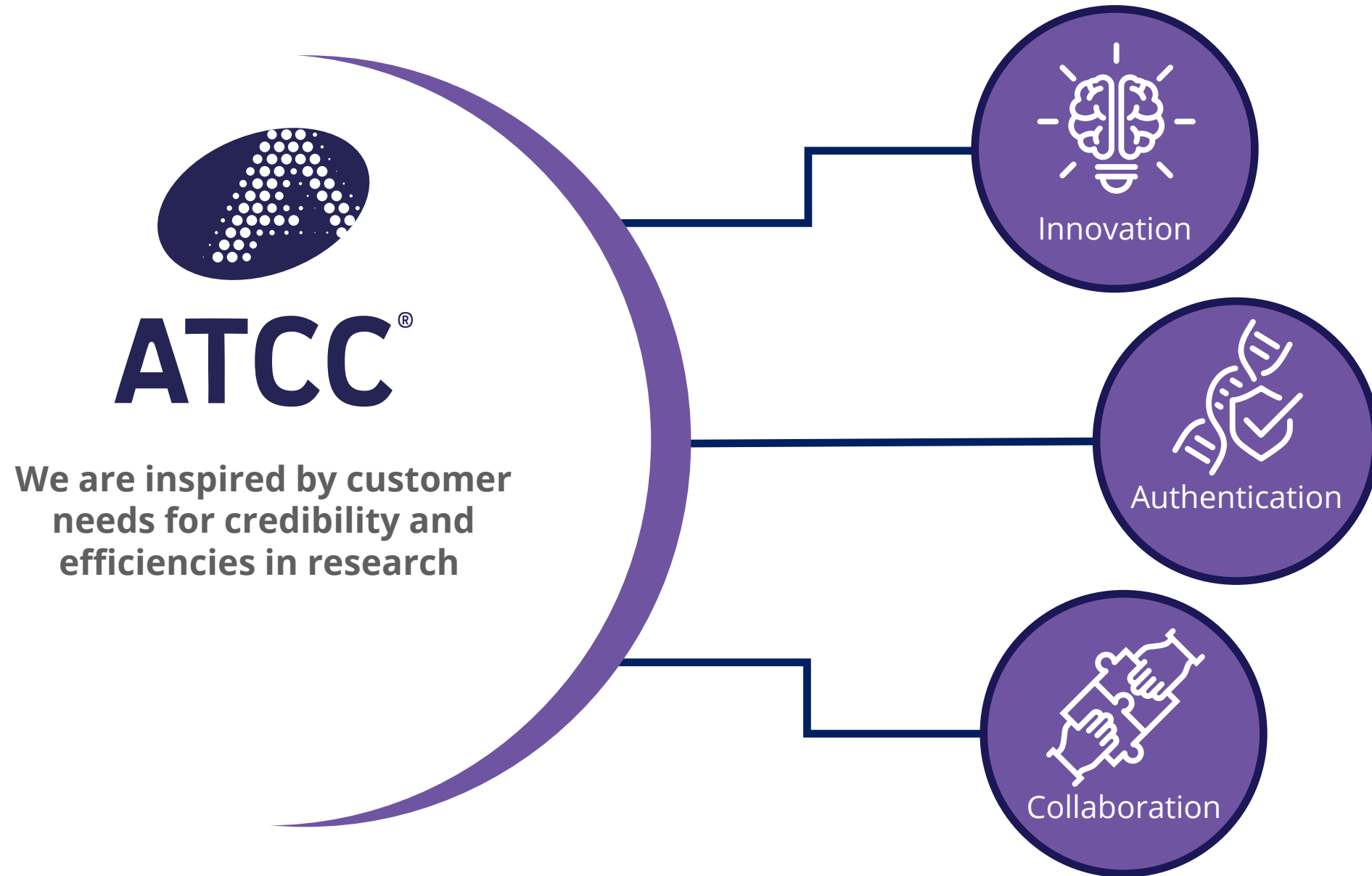

Microbial QC testing applications

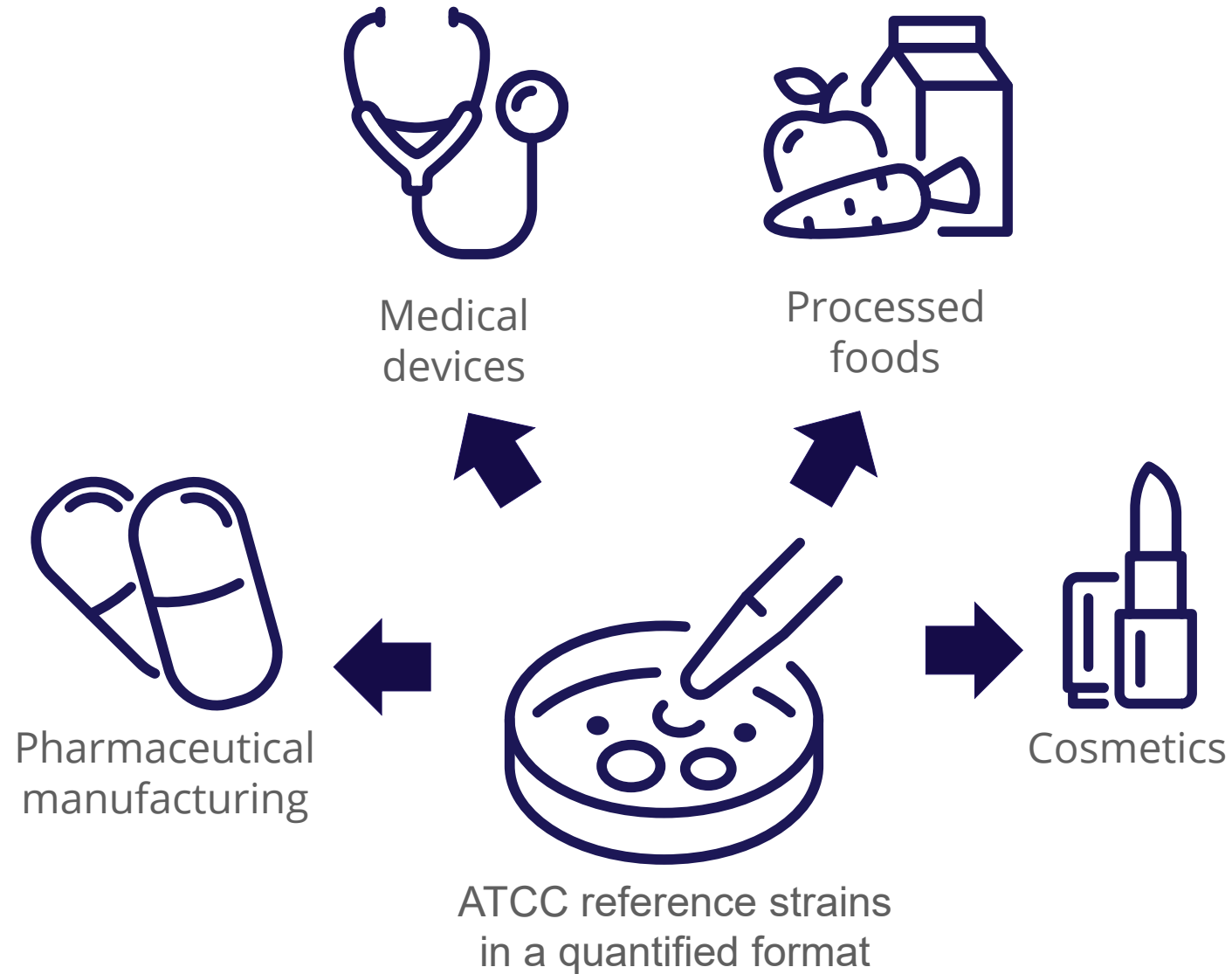

Pharmacopeias govern microbial QC testing

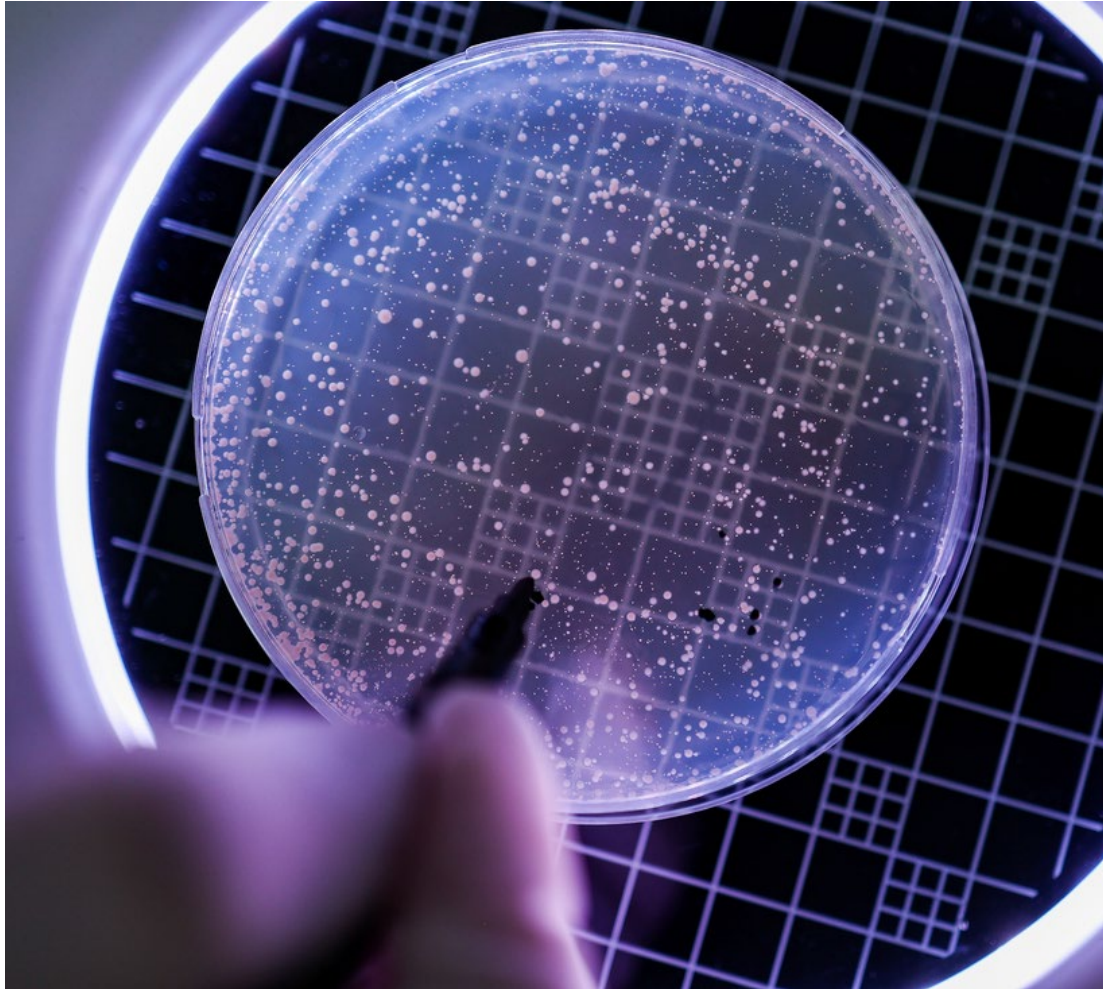


- Define standardized methods for microbial testing in pharmaceuticals and other regulated industries
- Ensure product safety, consistency, and regulatory compliance
- Provide for validated assays and acceptance criteria for microbial limits and preservative efficacy
- Used globally to guide assay development, assay validation, and routine quality controls

Challenges when using microbial reference strains



The traditional microbe culture workflow

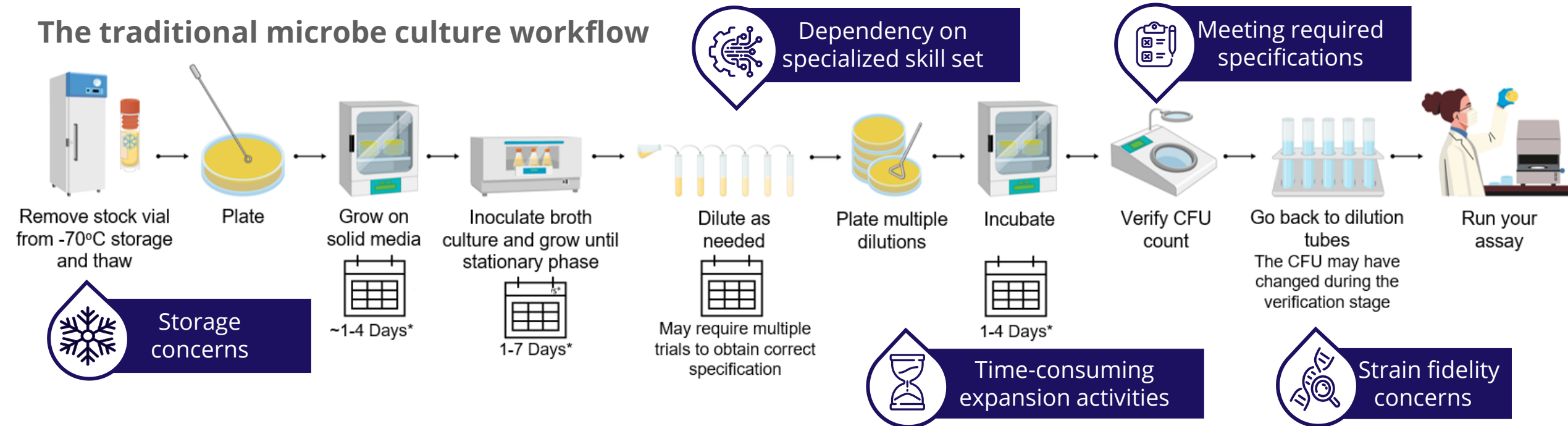

Challenges using single-use microbial reference strains



Features	Company A	Company B	Company C	Company D
Processing time ~1 minute			✓	✓
2-8°C refrigeration storage	✓	✓		
Manufactured under ISO 17034	✓	✓	✓	✓
Includes all strains for USP <51> (High CFU format)	✓		✓	✓
Includes all strains for USP <61> (Low CFU format)	✓	✓	✓	✓
Product kit includes rehydration buffer	✓	✓	✓	
Sourced from ATCC	✓	✓		
Provided as Passage 0 strain				

Challenges using single-use microbial reference strains



Features	MicroQuant™	Company A	Company B	Company C	Company D
Processing time ~1 minute	✓			✓	✓
2-8°C refrigeration storage	✓	✓	✓		
Manufactured under ISO 17034	✓	✓	✓	✓	✓
Includes all strains for USP <51> (High CFU format)	✓	✓		✓	✓
Includes all strains for USP <61> (Low CFU format)	✓	✓	✓	✓	✓
Product kit includes rehydration buffer	✓	✓	✓	✓	
Sourced from ATCC	✓	✓	✓		
Provided as Passage 0 strain	✓				




MicroQuant™
by **ATCC®**

Meeting the challenge through cryobiology




The traditional microbe culture workflow

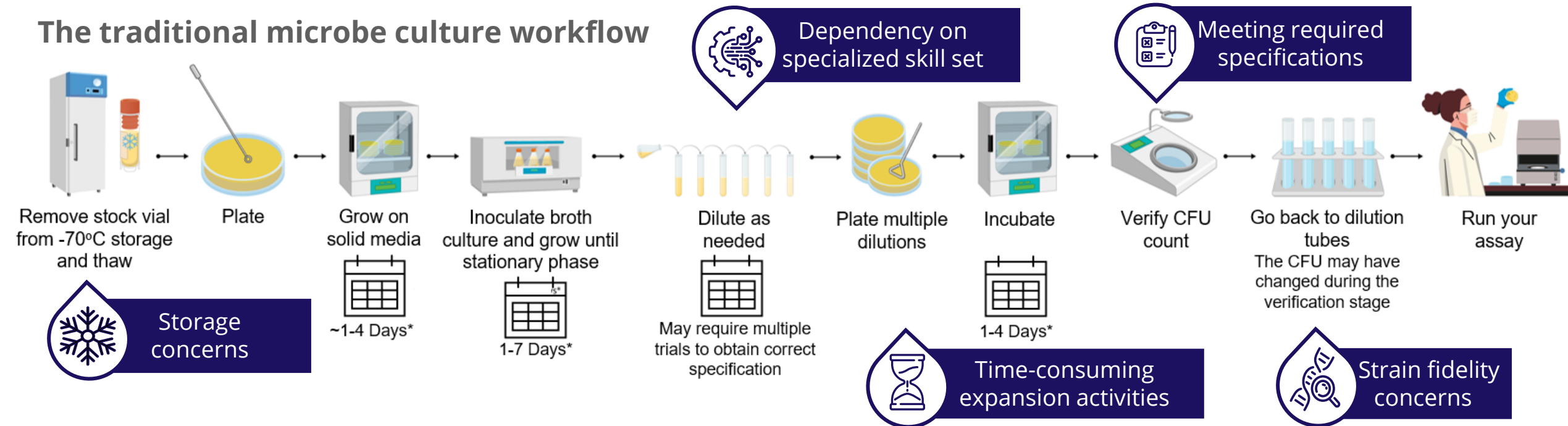


Meeting the challenge through cryobiology




The MicroQuant™ workflow

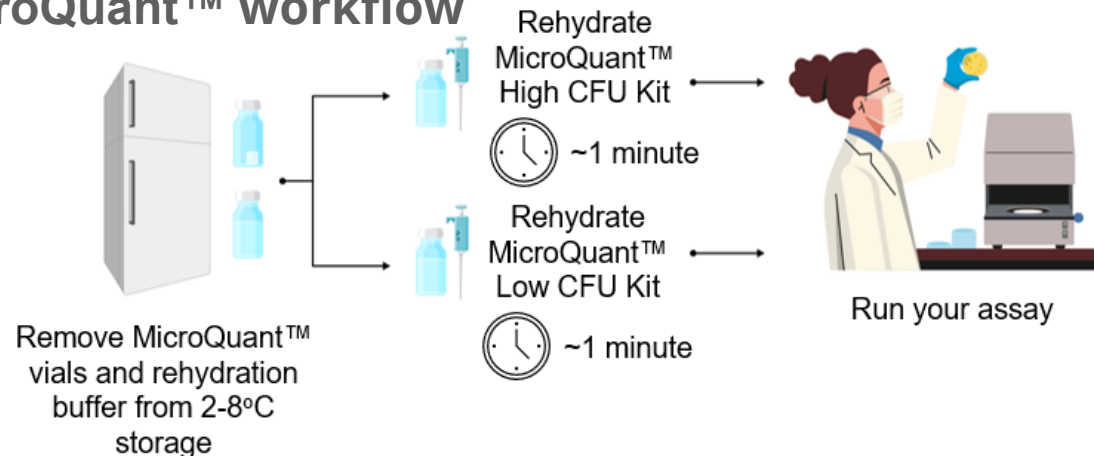

Our innovative cryopreservation technology enables you to streamline your microbial QC workflow from days to minutes.

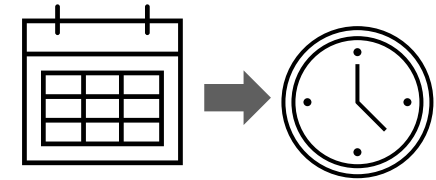

The traditional microbe culture workflow

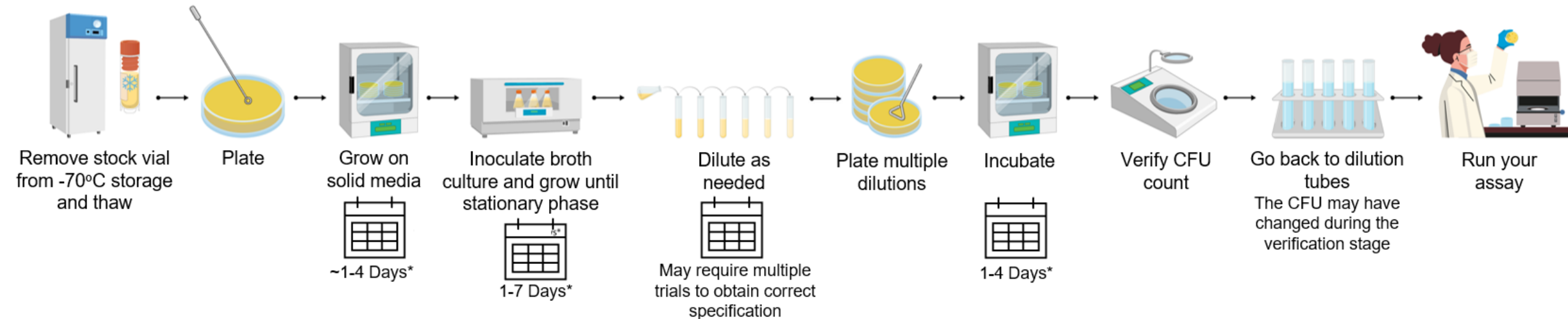



A ready-to-use solution to help you streamline microbial QC testing

- Precisely quantitated in high-titer and low-titer pelleted formats
- Single-use format using an innovative, proprietary preservation technology
- Rapid, uniform rehydration in less than a minute
- Easy to store and ready to use anytime—no need to thaw
- Original ATCC materials manufactured under ISO 17034

Precision backed by authentication




- **Phenotypic analysis** – Cell morphology, colony description, viability, and purity
- **Proteomic analysis** – MALDI-TOF MS
- **Functional testing** – Antimicrobial susceptibility testing, serotyping, virulence detection
- **Genotypic analysis** – Sequencing conserved regions of the genome and whole-genome sequencing
 - We have published the corresponding genomes for 5,500 ATCC microbes on our **ATCC® Genome Portal** (genomes.atcc.org)

Researchers look to ATCC for a wide range of authentication resources to safeguard reproducibility and meet requirements for funding, publication, and quality control.


MicroQuant™

by ATCC®

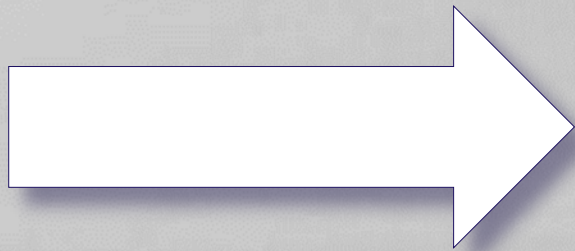

Maximize your resources

MicroQuant™ is stable at 2-8°C, making it convenient to store and ready to use anytime

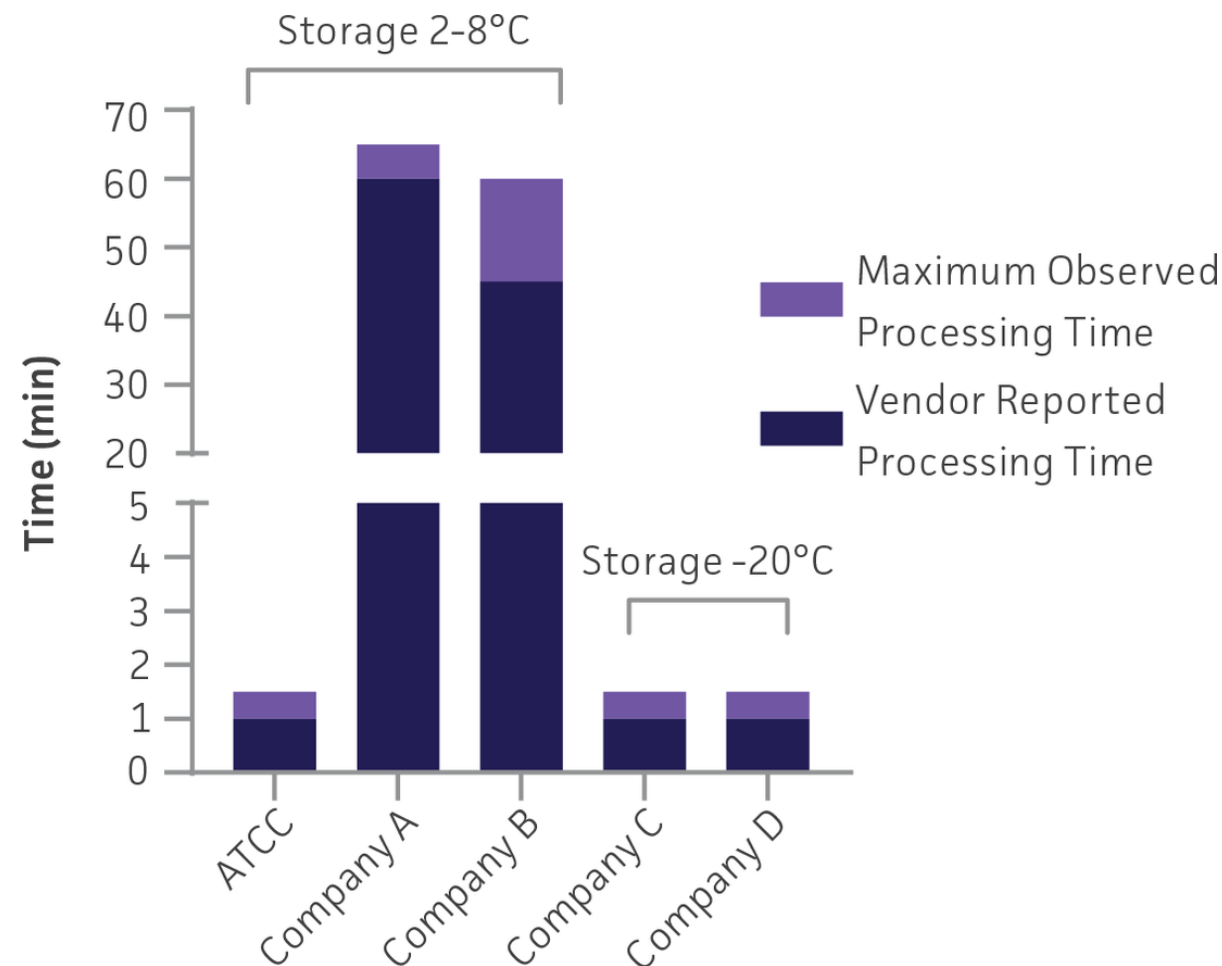


MicroQuant™ is stable up to 1 year at 2-8°C


Unlock flexibility with platform-agnostic solutions

MicroQuant™ is provided as a single-use quantitated pellet that enables fast assay set up with minimal handling

Low CFU Items
(100 to 1,000 CFU per pellet)

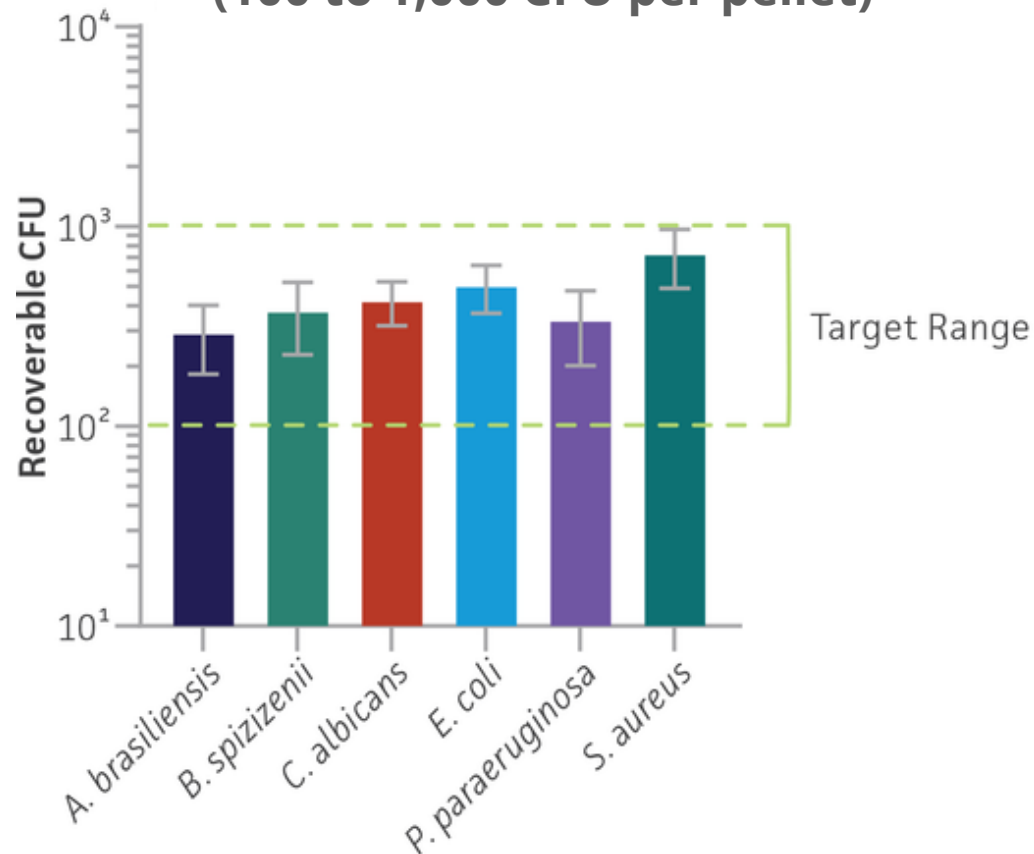







High CFU Items
(10^7 to 10^8 CFU per pellet)

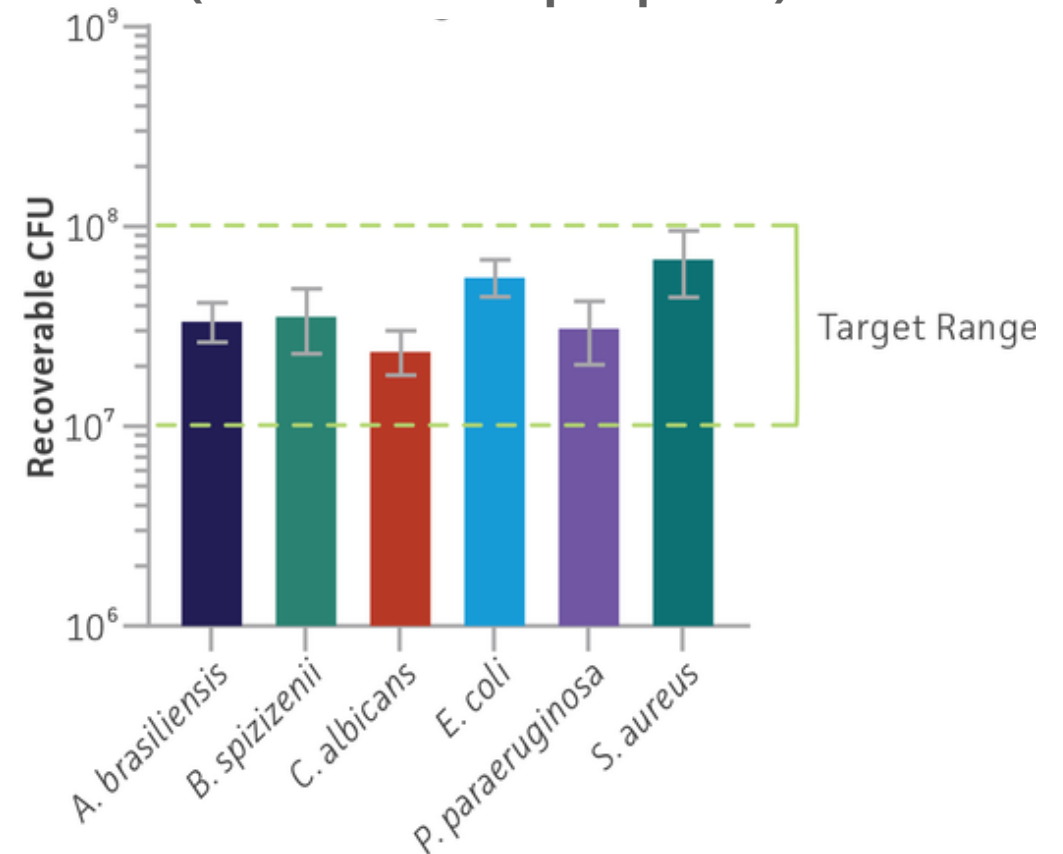

Rely on consistent, precise quantitation

MicroQuant™ - Precise and accurate low CFU pellets for bioburden testing







"Low CFU" means
100 to 1,000
CFU/pellet

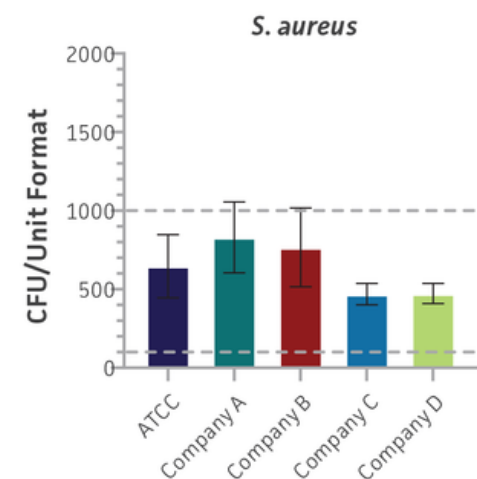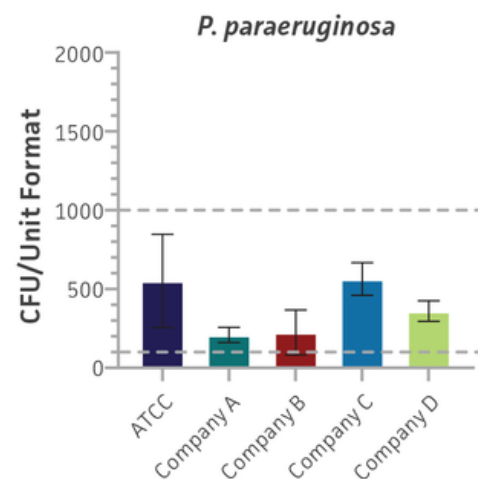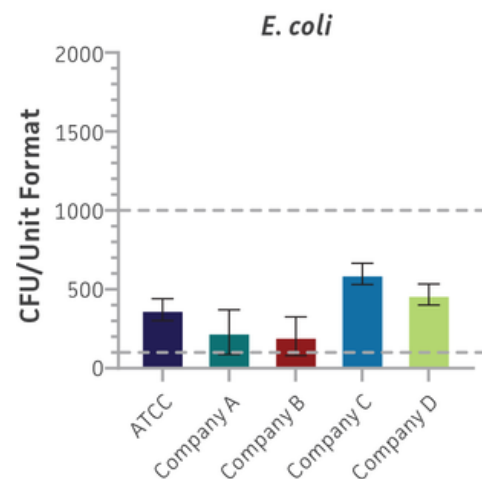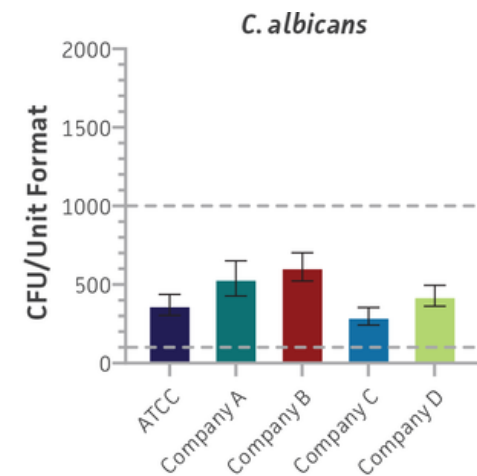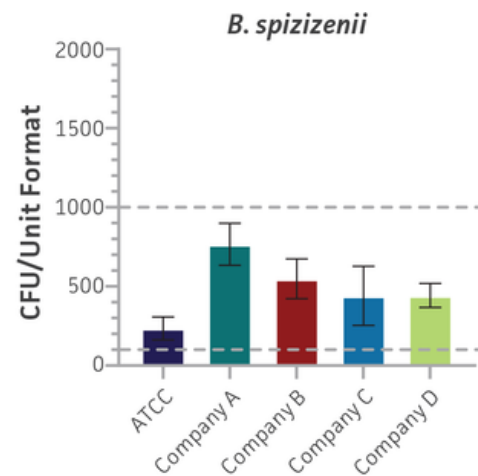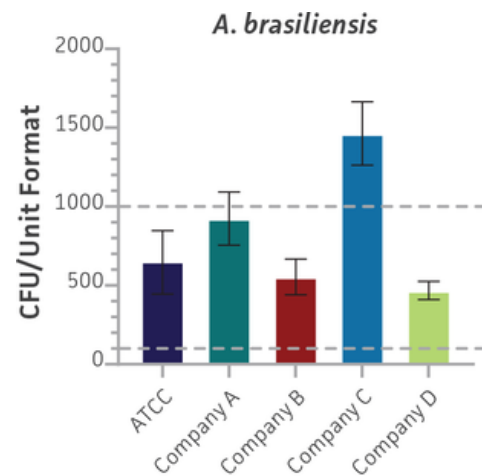

Rely on consistent, precise quantitation







MicroQuant™ - Precise, high CFU pellets for testing under USP <51> guidelines

"High CFU"
means
 10^7 to 10^8
CFU/pellet

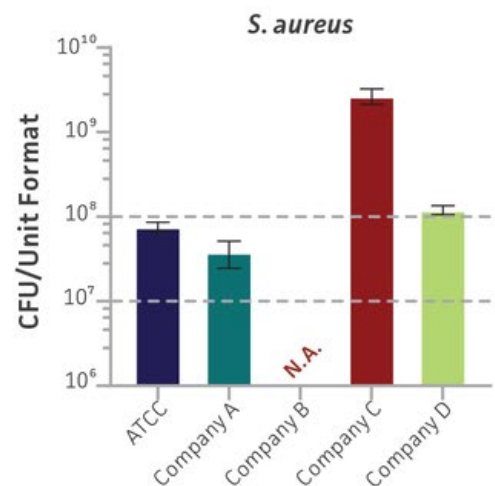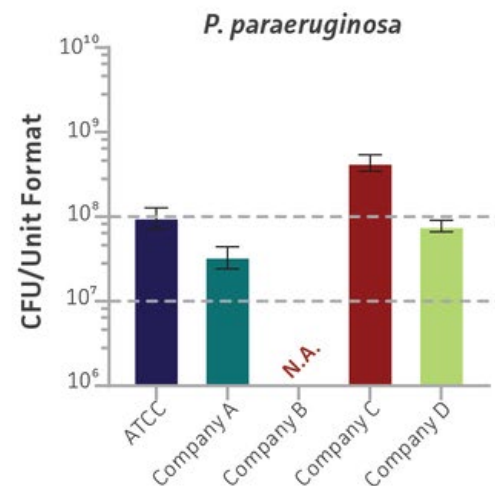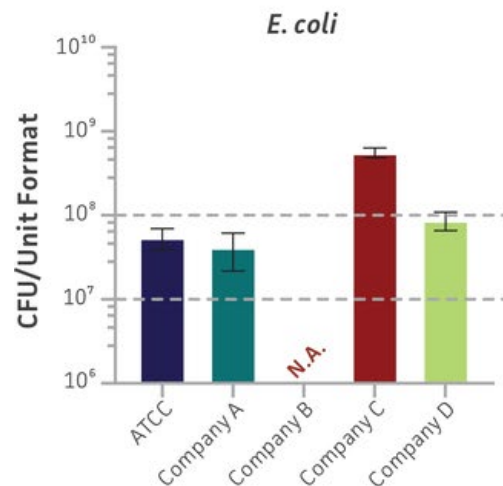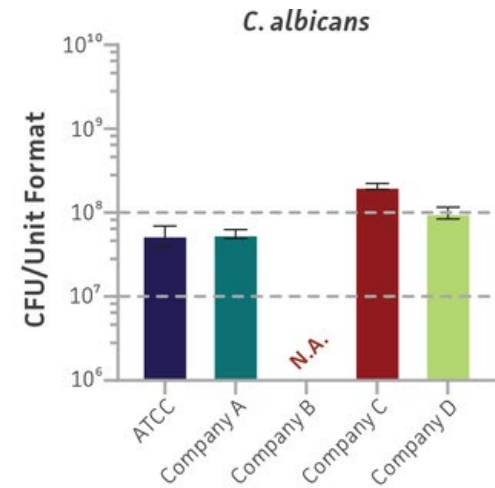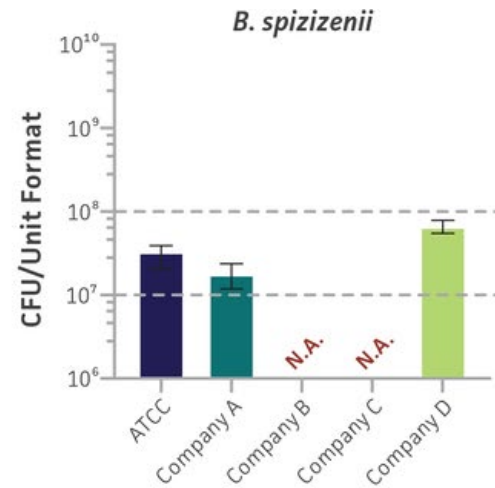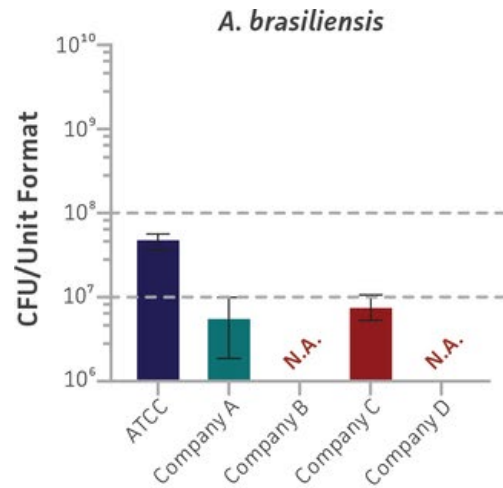







Short-term stability of MicroQuant™

Low titer

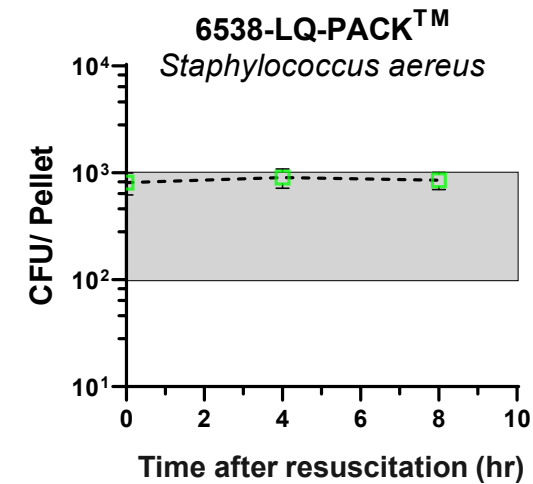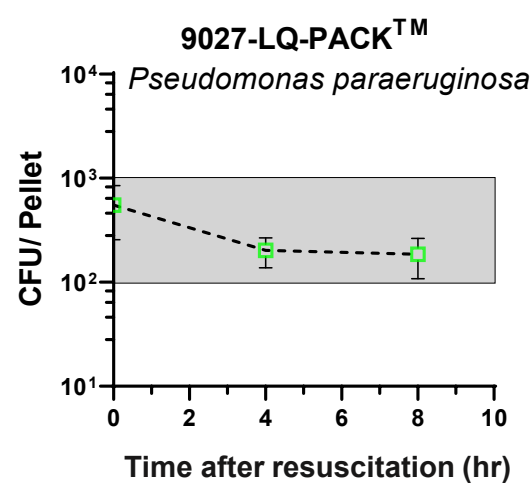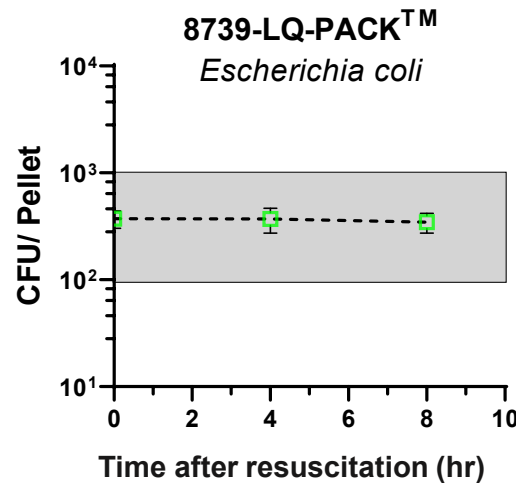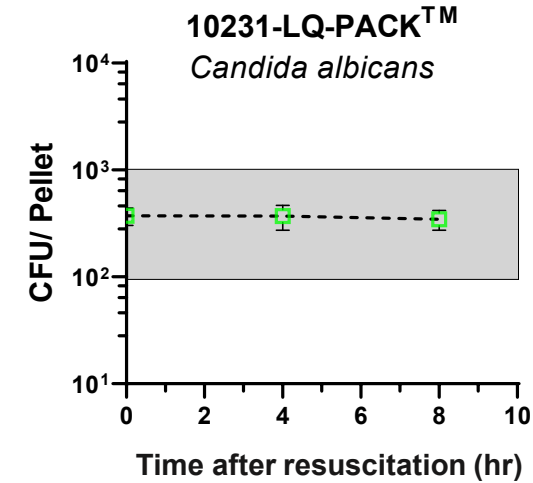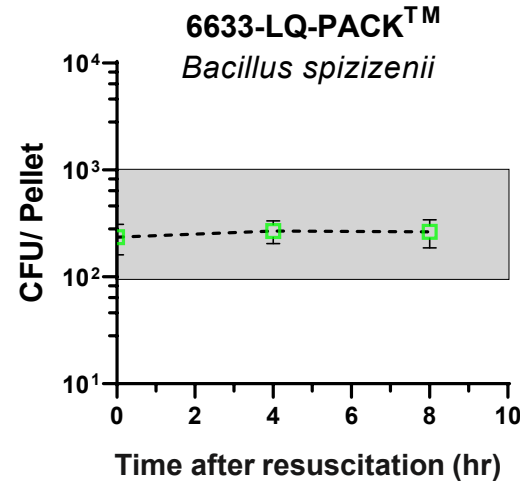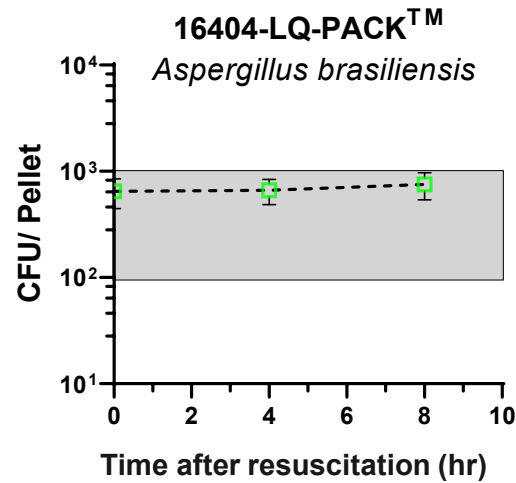







Short-term stability of MicroQuant™

High titer

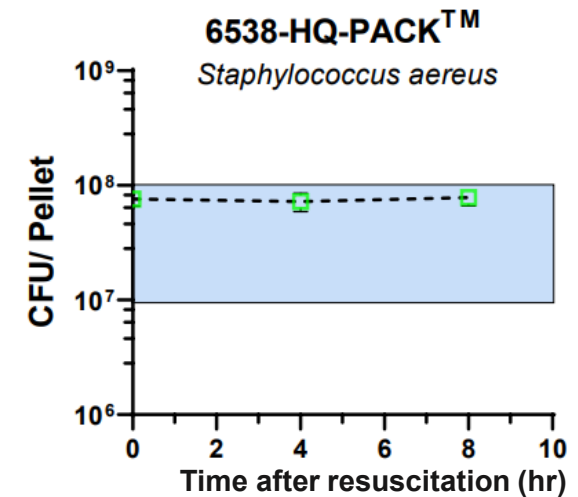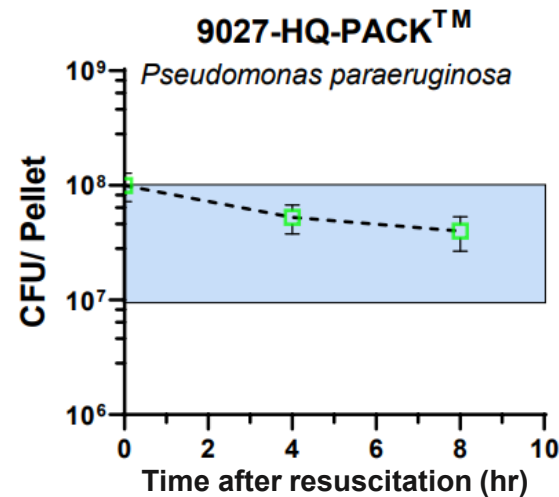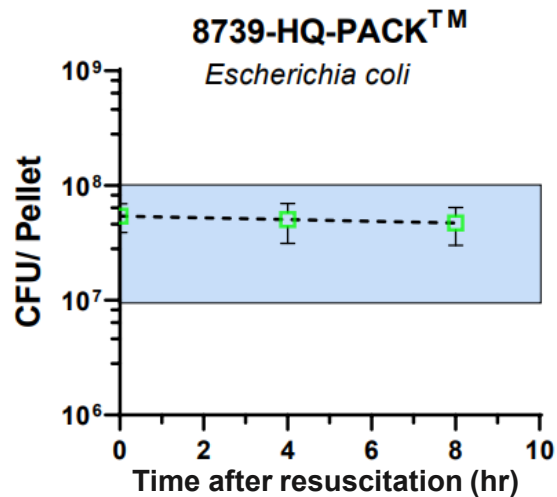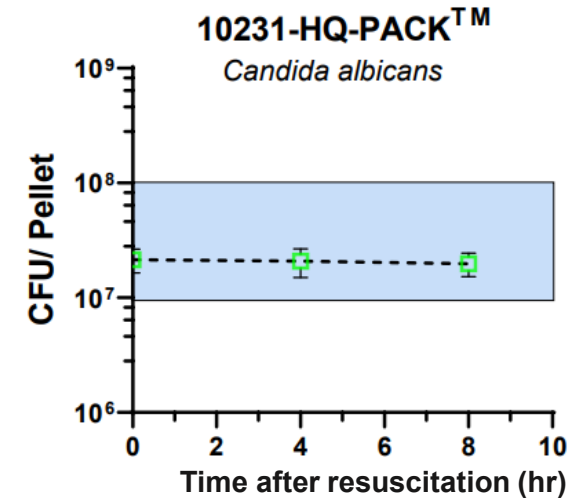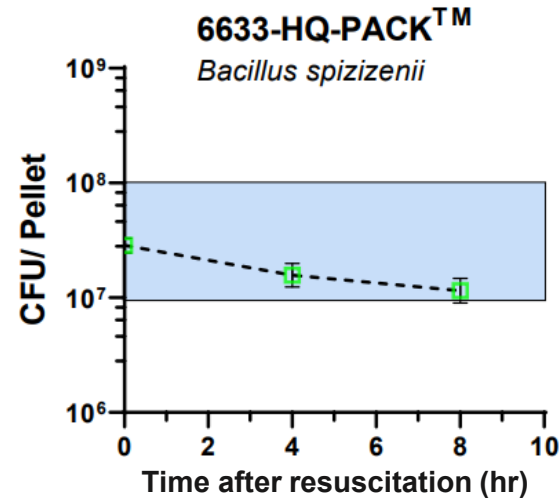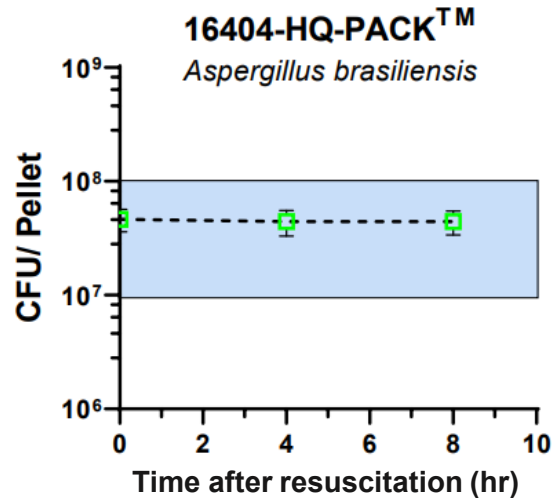


Transportation stability of MicroQuant™

Low titer

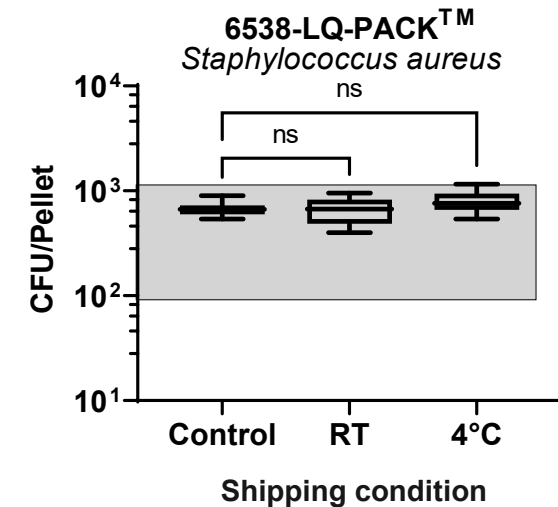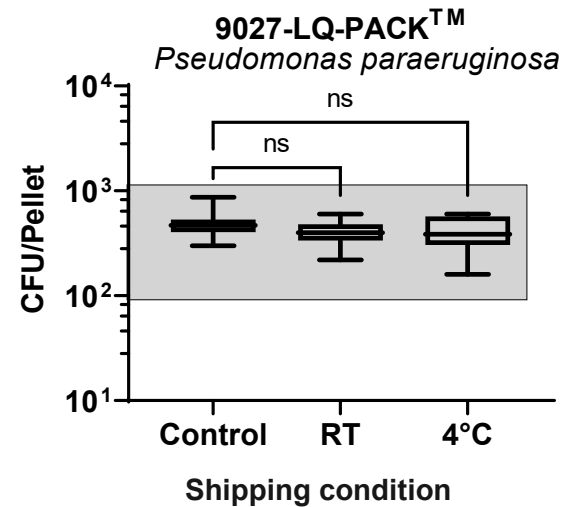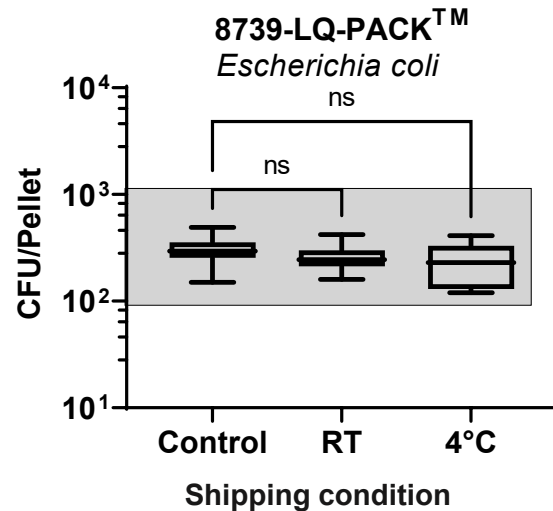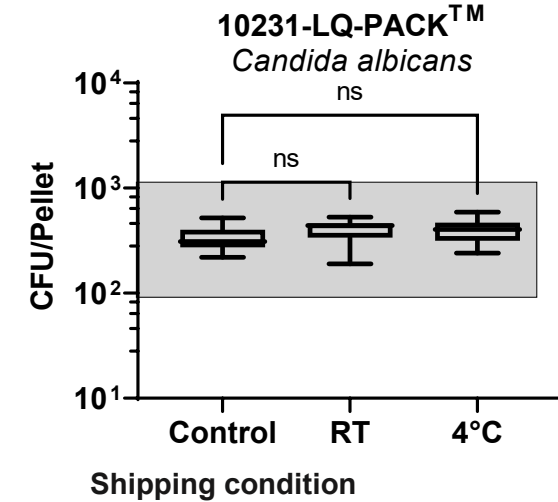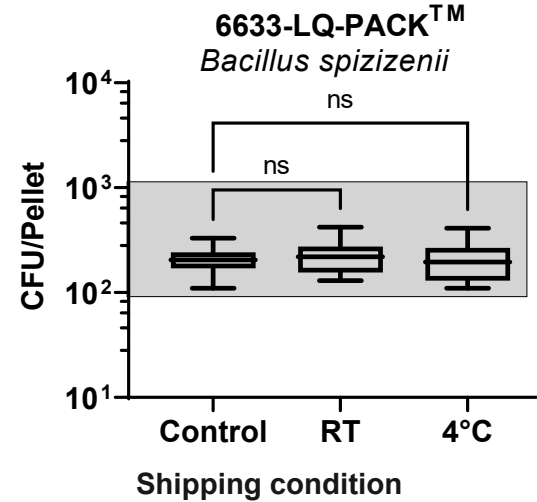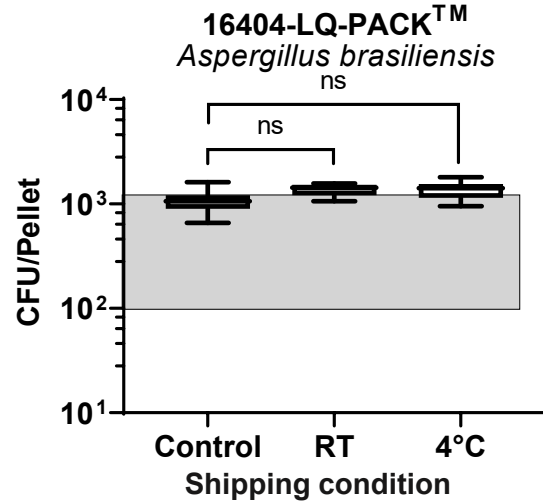


Transportation stability of MicroQuant™


High titer

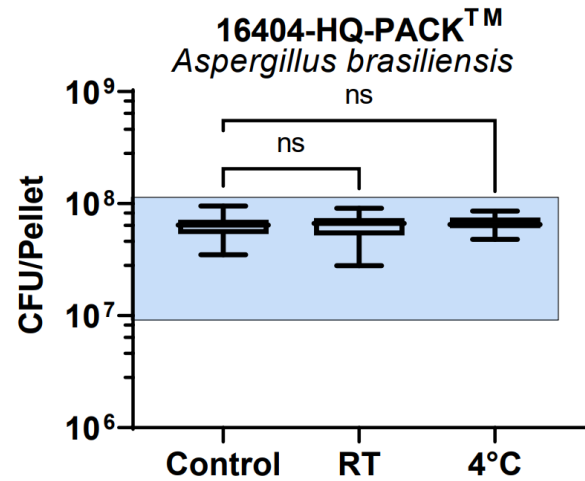


Shipping condition

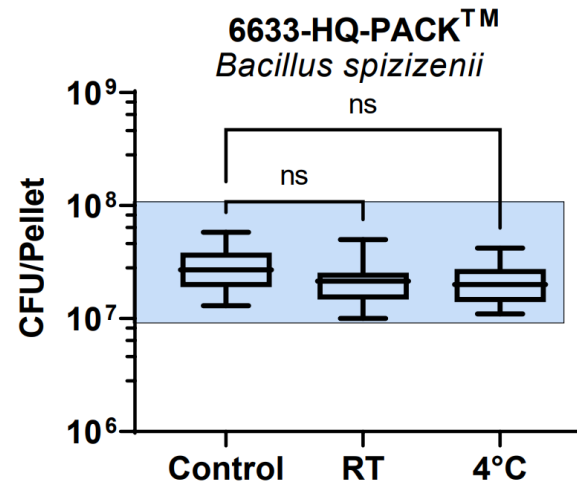


Shipping condition

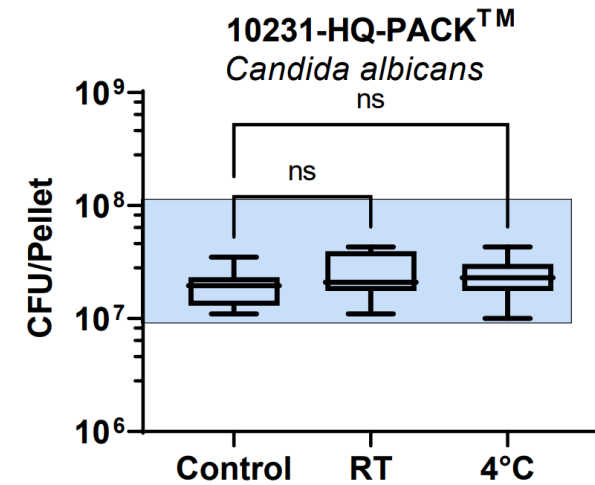


Shipping condition

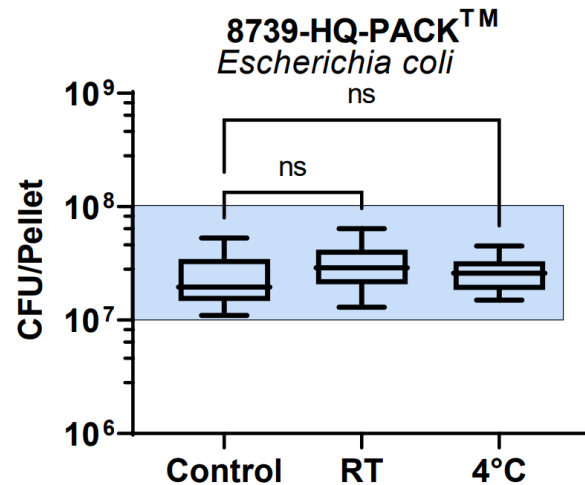

Shipping condition

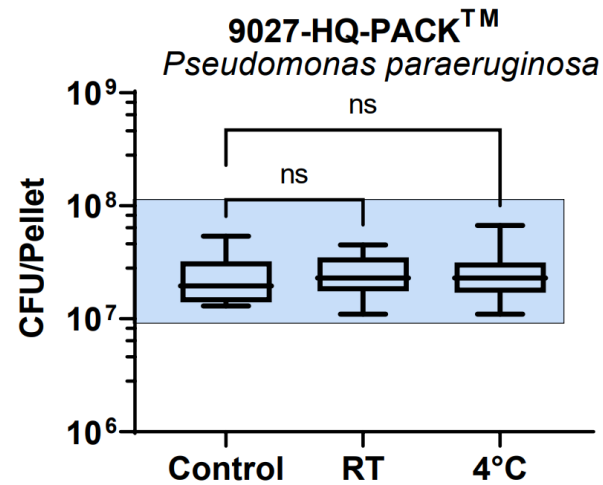

Shipping condition

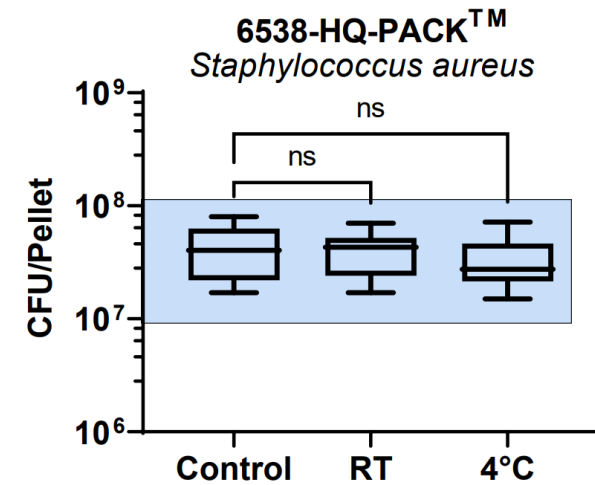

Shipping condition

How MicroQuant™ compares to similar products



Features	MicroQuant™	Company A	Company B	Company C	Company D
Processing time ~1 minute	✓			✓	✓
2-8°C refrigeration storage	✓	✓	✓		
Manufactured under ISO 17034	✓	✓	✓	✓	✓
Includes all strains for USP <51> (High CFU format)	✓	✓*		✓*	✓*
Includes all strains for USP <61> (Low CFU format)	✓	✓	✓	✓	✓
Product kit includes rehydration buffer	✓	✓	✓	✓	
Sourced from ATCC	Passage 0	✓	✓		

*Some strains were not available for purchase at the time of testing

End-user benefits and feedback on using MicroQuant™

MicroQuant™ saves time and cost

Versus preparing internal controls



86%

of labor time is saved when using MicroQuant™ versus preparing internal controls manually

83%

cost reduction is achieved using MicroQuant™ versus using internally prepared controls



External end-user results for quantitation

MicroQuant™ is a top performer compared to other single-use controls

End-user verification of reported CFU/mL for single-use control per Instructions for Use

Single-use control	Strains used in high titer (USP <51>)			Strains used in low titer (USP <61>)		
	<i>C. albicans</i>	<i>S. aureus</i>	<i>E. coli</i>	<i>C. albicans</i>	<i>S. aureus</i>	<i>B. spizizenii</i>
MicroQuant™ by ATCC	✓	✓	✓	✓	✓	✓
Company A	✓	✓	✓	✓	✓	✓
Company B	✓	✓	✓	X	X	X
Company C	Not tested	Not tested	Not tested	✓	Not tested	✓
Company D	✓	X	X	✓	✓	X

✓ = met criteria for product's instruction of use
 X = didn't meet criteria for product's instruction of use

MicroQuant™ had the "most accurate CFU results"

External end-user results for ease of dissolution

MicroQuant™ is the top performer compared to other single-use controls



Ease of dissolution scored on a 1 to 5 rating

Single-use control	<i>C. albicans</i> / <i>S. aureus</i> / <i>E.coli</i>
MicroQuant™ by ATCC	★★★★★
Company A	★★☆☆☆
Company B	★★★★☆
Company C	★★☆☆☆
Company D	★★☆☆☆

Score: 1 to 5 system

- 5: Fully soluble within seconds at room temperature.
- 4: Dissolves quickly (<1 minute) with minimal effort.
- 3: Dissolves with moderate effort (10 minutes, stirring).
- 2: Requires extended processing or specific steps to dissolve.
- 1: Insoluble or requires elevated temperatures/incubation.

MicroQuant™
had the
“fastest
rehydration”

External end-user results for ease of preparation

MicroQuant™ is the top performer compared to other single-use controls



Ease of preparation scored on a 1 to 5 rating

Single-use control	Observation rating	Justification
MicroQuant™ by ATCC	★★★★★	Simple and intuitive preparation. No clumping or precipitation. Buffer included.
Company A	★★★☆☆	Moderate preparation with 1 hour of warming . Additional buffer needed for some formats.
Company B	★★★☆☆	Simple process but physical risks noted (pellet may miss vial, tipping risk).
Company C	★★★★☆	Simple preparation. No consumables needed. Incubator required.
Company D	★★★★☆	Simple preparation. No clumping or precipitation noted. Additional buffer needed for high CFU quantitation.

MicroQuant™
had the most
“intuitive prep”

What MicroQuant™ products are available?

Explore the MicroQuant™ portfolio



ATCC® No.	Description
10231-LQ-PACK™	MicroQuant™ <i>Candida albicans</i> , low CFU
10231-HQ-PACK™	MicroQuant™ <i>Candida albicans</i> , high CFU
9027-LQ-PACK™	MicroQuant™ <i>Pseudomonas paraeruginosa</i> , low CFU
9027-HQ-PACK™	MicroQuant™ <i>Pseudomonas paraeruginosa</i> , high CFU
6538-LQ-PACK™	MicroQuant™ <i>Staphylococcus aureus</i> subsp. <i>aureus</i> , low CFU
6538-HQ-PACK™	MicroQuant™ <i>Staphylococcus aureus</i> subsp. <i>aureus</i> , high CFU
16404-LQ-PACK™	MicroQuant™ <i>Aspergillus brasiliensis</i> , low CFU
16404-HQ-PACK™	MicroQuant™ <i>Aspergillus brasiliensis</i> , high CFU
8739-LQ-PACK™	MicroQuant™ <i>Escherichia coli</i> , low CFU
8739-HQ-PACK™	MicroQuant™ <i>Escherichia coli</i> , high CFU
6633-LQ-PACK™	MicroQuant™ <i>Bacillus spizizenii</i> , low CFU
6633-HQ-PACK™	MicroQuant™ <i>Bacillus spizizenii</i> , high CFU

Low CFU = 100 to 1,000 CFU/pellet

High CFU = 10^7 to 10^8 CFU/pellet



Explore the MicroQuant™ portfolio



ATCC® No.	Product Name	Description
MQ-51™	MicroQuant™ Antimicrobial Effectiveness Panel	<p>A panel comprising the five microbial challenge organisms cited in USP <51>.</p> <p>Each component is provided as a pack containing 5 vials of cryopreserved pellets (10^7 to 10^8 CFU per pellet) and 5 vials of rehydration buffer.</p>
MQ-61™	MicroQuant™ Microbial Examination of Nonsterile Products Panel	<p>A panel comprising the five microbial challenge organisms cited in USP <61>.</p> <p>Each component is provided as a pack containing 5 vials of cryopreserved pellets (100 to 1,000 CFU per pellet, supporting 10 assays per pellet) and 5 vials of rehydration buffer.</p>



MicroQuant™ applications

Control strains supporting USP general chapter requirements



High-quantitation (HQ)
Low-quantitation (LQ)

Species		USP <51>	USP <61>	USP <60>	USP <62>	USP <71>	USP <72>
Available Now	<i>Aspergillus brasiliensis</i> WLRI 034(120)	HQ	LQ			LQ	LQ
	<i>Bacillus spizizenii</i> NRS 231		LQ			LQ	LQ
	<i>Candida albicans</i> 3147	HQ	LQ		LQ	LQ	LQ
	<i>Escherichia coli</i> Crooks	HQ			LQ		
	<i>Pseudomonas paraeruginosa</i> R. Hugh 813	HQ	LQ	LQ	LQ	LQ	LQ
	<i>Staphylococcus aureus</i> FDA 209	HQ	LQ	LQ	LQ	LQ	LQ
Available in July	<i>Burkholderia cepacia</i> UCB 717			LQ			
	<i>Burkholderia cenocepacia</i> LMG 16656			LQ			
	<i>Burkholderia multivorans</i> LMG 13010			LQ			
Available by the end of 2025	<i>Clostridium sporogenes</i> L.S. McClung 2006				LQ	LQ	LQ
	<i>Clostridium sporogenes</i> SR 5				LQ	LQ	LQ
	<i>Salmonella enterica</i> CDC 6516-60				LQ		

Summary



- ATCC® MicroQuant™ is an innovative product suite designed to streamline microbial quality control testing.
- Precisely quantitated
 - High-titer (HQ; 10^7 to 10^8 CFU per pellet)
 - Low-titer (LQ; 100 to 1,000 CFU per pellet)
- Single-use format for fast assay setup and minimal handling.
- Developed from traceable, original source materials provided at **Passage zero**.
- Immediate rehydration at room temperature with an 8-hour usability window.
- Stable storage at 2-8°C.
- Manufactured under an ISO 17034–accredited process.





ATCC[®]

CREDIBLE LEADS TO INCREDIBLE

Thank You