

**ATCC Medium 274: Tryptone Broth/ Agar/ Stabs, 1%**

**Broth**

Tryptone..... 10.0 g  
DI water..... 1000.0 mL

Mix well. Dispense as required and autoclave at 121°C for 15 minutes.

**Agar**

Tryptone..... 10.0 g  
DI water..... 1000.0 mL  
Agar..... 15.0 g

Mix well. Autoclave at 121°C for 15 minutes and dispense as required.

**Stabs**

Tryptone..... 10.0 g  
DI water..... 1000.0 mL  
Agar (0.5%)..... 5.0 g

Mix well and boil to dissolve agar. Dispense as required and autoclave at 121°C for 15 minutes.