

**ATCC medium: 1334 TPL medium**

Potato (see below).....	200.0 g
Yeast extract.....	30.0 g
Liver (see below).....	25.0 g
Beef extract.....	5.0 g
Thioglycollate medium (dehydrated).....	10.0 g
Glucose.....	5.0 g
Glycerol.....	15.0 g
Agar.....	15.0 g
Distilled water to.....	1.0 L

Adjust to pH 7.0.

Gently boil sliced potatoes in 500 ml of the water (or sliced liver in 150 ml of the water) for 30 minutes and remove solids by filtration through cloth.

Autoclave at 121C for 15 minutes.