



## NDM-1 STRAINS

In the last decade, confirmed NDM-1 strains have been attributed to numerous nosocomial infections worldwide. Regrettably, the isolation and identification of NDM-1 strains is time consuming and often not routinely available in clinical laboratories, further contributing to escalating mortality rates.

ATCC understands the danger and growing concern behind the spread of multidrug-resistance strains. To aid in the prevention and treatment of NDM-1-associated infections, ATCC has acquired and authenticated multidrug-resistant prokaryotic isolates known to produce NDM-1.

Each clinical strain has undergone biochemical testing, extensive antibiotic profiling, and genomic analysis to verify:

- Species authentication
- NDM-1 production
- Multidrug-resistance

The table provided below lists the NDM-1 strains available from ATCC.

### ATCC® NDM-1 Strains

ATCC® No.	Strain Description	Strain Designation	Presence of Select Virulence Genes
<a href="#">BAA-2146™</a>	<i>Klebsiella pneumoniae</i>	1000527, 7561	<i>bla</i> <sub>NDM</sub> +/ <i>bla</i> <sub>KPC</sub> -
<a href="#">BAA-2452™</a>	<i>Escherichia coli</i>	NDM-1	<i>bla</i> <sub>NDM</sub> +/ <i>bla</i> <sub>KPC</sub> -
<a href="#">BAA-2468™</a>	<i>Enterobacter cloacae</i>	1000654	<i>bla</i> <sub>NDM</sub> +/ <i>bla</i> <sub>KPC</sub> -
<a href="#">BAA-2469™</a>	<i>Escherichia coli</i>	1001728	<i>bla</i> <sub>NDM</sub> +/ <i>bla</i> <sub>KPC</sub> -
<a href="#">BAA-2470™</a>	<i>Klebsiella pneumoniae</i> subsp. <i>pneumoniae</i>	1002565	<i>bla</i> <sub>NDM</sub> +/ <i>bla</i> <sub>KPC</sub> -
<a href="#">BAA-2471™</a>	<i>Escherichia coli</i>	1100101	<i>bla</i> <sub>NDM</sub> +/ <i>bla</i> <sub>KPC</sub> -
<a href="#">BAA-2472™</a>	<i>Klebsiella pneumoniae</i> subsp. <i>pneumoniae</i>	1100975	<i>bla</i> <sub>NDM</sub> +/ <i>bla</i> <sub>KPC</sub> -
<a href="#">BAA-2473™</a>	<i>Klebsiella pneumoniae</i>	1100770	<i>bla</i> <sub>NDM</sub> +/ <i>bla</i> <sub>KPC</sub> -

### ATCC® NDM-1 Panel

The strains are available individually, or grouped together in the ATCC® NDM-1 Strains Panel (ATCC® MP-18™) to allow for convenience in ordering and bulk discount pricing. The strains are ideal for the development, verification, and evaluation of:

- Rapid detection methods
- Innovative therapeutic techniques
- Novel antibiotics
- Updated sterility protocols

ATCC hopes that with your diligent research and our authenticated strains, we can stop the spread of these superbugs together!

### Background on NDM-1

Gram-negative prokaryotes expressing New Delhi metallo-β-lactamase (NDM-1), an enzyme that confers resistance to beta-lactam antibiotics including last-resort carbapenem antibiotics, are frequently multi-drug resistant. NDM-1 is encoded by the plasmid-associated gene, *bla*<sub>NDM</sub>. Horizontal transfer of this mobile genetic element has allowed NDM-1 to be passed between Enterobacteriaceae genera commonly found in the human microbiome, including *Escherichia coli*, *Enterobacter cloacae*, and *Klebsiella pneumoniae*. The spread of these multidrug-resistant strains can be further enhanced by inadvertent transmission between immunocompromised patients in critical care facilities or through the transfer of infected patients between healthcare facilities.

### ATCC® NDM-1 Panel (MP-18™)

Take advantage of convenient ordering and bulk discounts associated with purchasing an ATCC panel. The ATCC® NDM-1 Panel (MP-18™) is comprised of the following strains:

- ATCC® [BAA-2468™](#)
- ATCC® [BAA-2469™](#)
- ATCC® [BAA-2470™](#)
- ATCC® [BAA-2471™](#)
- ATCC® [BAA-2472™](#)
- ATCC® [BAA-2473™](#)

Enhance your research today with ATCC® [MP-18™](#).

ATCC® NDM-1 Strains – Antibiotic Profiles

		<a href="#">BAA-2146™</a>	<a href="#">BAA-2452™</a>	<a href="#">BAA-2468™</a>	<a href="#">BAA-2469™</a>	<a href="#">BAA-2470™</a>	<a href="#">BAA-2471™</a>	<a href="#">BAA-2472™</a>	<a href="#">BAA-2473™</a>
Penicillins	Amoxicillin/ Clavulanic Acid	R	R	R	R	R	R	R	R
	Ticarcillin	R	R	R	R	R	R	R	R
	Ticarcillin/ Clavulanic Acid	R	Not Tested	R	R	R	R	R	R
	Piperacillin	R	R	R	R	R	R	R	R
	Ampicillin	R	R	R	R	R	R	R	R
	Ampicillin/ Sulbactam	R	R	R	R	R	R	R	R
Cephalosporins	Cefalotin	R	R	R	R	R	R	R	R
	Cefuroxime	R	R	R	R	R	R	R	R
	Cefuroxime Axetil	R	R	R	R	R	R	R	R
	Cefotetan	R	R	R	R	R	R	R	R
	Cefpodoxime	R	R	R	R	R	R	R	R
	Cefotaxime	R	R	R	R	R	R	R	R
	Ceftizoxime	R	R	R	R	R	R	R	R
	Cefazolin	R	R	R	R	R	R	R	R
	Cefoxitin	R	R	R	R	R	R	R	R
	Ceftazidime	R	R	R	R	R	R	R	R
	Ceftriaxone	R	R	R	R	R	R	R	R
Cefepime	R	R	R	R	R	R	R	R	
Carbapenems	Doripenem	R	Not Tested	R	R	R	R	R	R
	Meropenem	R	R	R	R	R	R	R	R
	Ertapenem	R	R	R	R	R	R	R	R
	Imipenem	R	R	R	R	R	R	S	R
Quinolones	Nalidixic Acid	R	S	R	R	R	R	R	R
	Moxifloxacin	R	S	R	R	R	R	R	R
	Norfloxacin	R	S	R	R	R	R	R	R
	Ciprofloxacin	R	S	R	R	R	R	R	R
	Levofloxacin	R	S	R	R	R	R	R	R
Amino-glycosides	Amikacin	R	R	R	R	R	I	R	R
	Gentamicin	R	R	R	R	R	R	R	R
	Tobramycin	R	R	R	R	R	R	R	R
Other	Tetracycline	R	S	R	R	R	R	R	R
	Tigecycline	R	S	R	S	S	S	R	R
	Nitrofurantoin	R	S	R	S	R	I	R	R
	Aztreonam	R	R	R	R	R	R	R	R
	Trimethoprim/ Sulfamethoxazole	R	S	R	R	R	R	R	R

R = Resistant, S = Susceptible, I = Intermediate susceptibility

**PHONE**  
 800.638.6597  
 703.365.2700

**EMAIL**  
 sales@atcc.org  
 tech@atcc.org

**WEB**  
 www.atcc.org



Learn more about ATCC Microbial Panels at [www.atcc.org/mp](http://www.atcc.org/mp)

10801 University Blvd.  
 Manassas, VA 20110

MB-1215-03

©2015 American Type Culture Collection. The ATCC trademark and trade name, and any other trademarks listed in this publication are trademarks owned by the American Type Culture Collection unless indicated otherwise. These products are for laboratory use only. Not for human or diagnostic use. ATCC products may not be resold, modified for resale, used to provide commercial services or to manufacture commercial products without prior ATCC written approval.